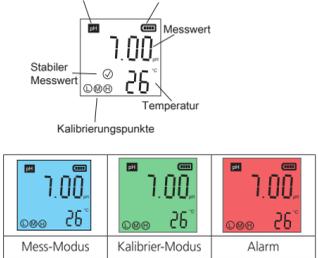
Display

Mess-Modus Kalibrier-Modus Alarm

Tastaturfunktion

Kurzer Tastendruck (1 Sek.) - Langer Tastendruck (3 Sek.)

Table with 3 columns: Action (Short/Long press), Mode (Mess/Aus/Calib), and Function (e.g., power on/off, backlight, settings, calibration).

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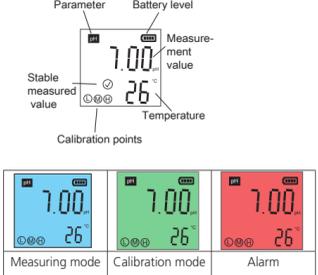
Display

Measuring mode Calibration mode Alarm

Keyboard function

Short keystroke (1 sec.) - Long keystroke (3 sec.)

Table with 3 columns: Action (Short/Long press), Mode (measuring/off/calib), and Function (power, backlight, settings, calibration).

Settings

In off mode, press 3 seconds: Opening Settings.


Navigate Confirm

Blue marked = factory setting

Table with 3 columns: Parameter (pH calibration series, alarm limit), Selection (USA, NIST, pH 0-14), and Meaning (description).

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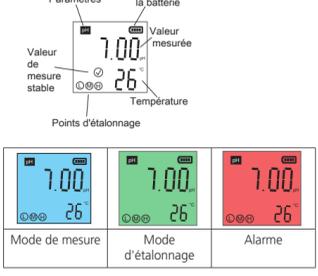Pression courte (1 sec.) - Pression longue (3 sec.)

Table with 3 columns: Action (Short/Long press), Mode (mesure/arrêt/calib), and Function (power, backlight, settings, calibration).

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
pH Kalibrierung

Table with 4 columns: Step number, Action (immersion, button press), and Description (1-point, 2-point, 3-point calibration).

Table with 3 columns: Parameter (Auto-Hold, Backlight, Temperature unit, Factory calibration), Value (On/Off, 1 min, °C/°F, Yes/No), and Description.

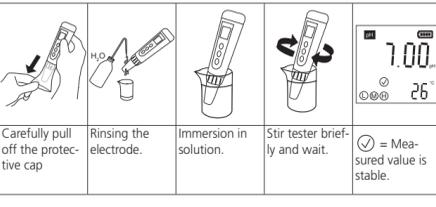pH measurement

Note: Press at "Hold" for new measurement recording.

pH Calibration


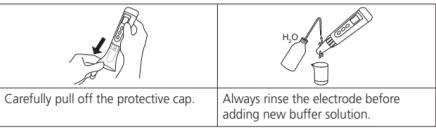
Table with 4 columns: Step number, Action (immersion, button press), and Description (1-point, 2-point, 3-point calibration).

After successful 1, 2 or 3 point calibration, the display shows:

Table showing display icons for 1-point, 2-point, and 3-point calibration.

Table with 3 columns: Parameter (Max. limit alarm, Auto-Hold, backlight, temperature unit, factory calibration), Value (pH 0-14, On/Off, 1 min, °C/°F, Yes/No), and Description.

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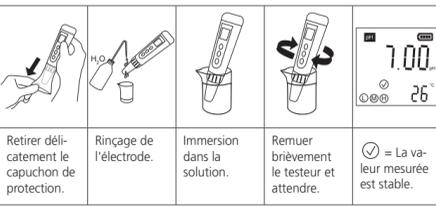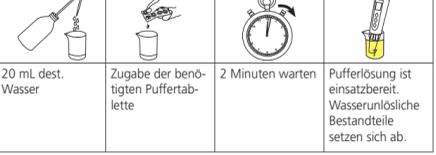
Austausch Batterien


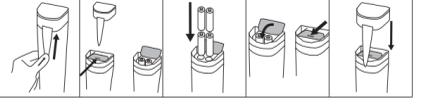
Table showing display icons for 3-point calibration and pH 4/7/10.

Note: Other pH calibration series possible according to settings.

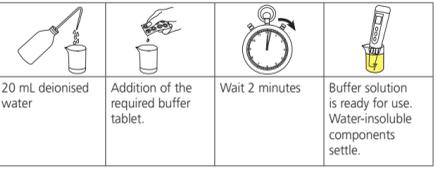Preparation of the buffer solution


Note: Use the buffer solution only once. The maximum deviation from the specified pH value is ± 0.05 pH.

SD 10 pH/T Specifications

Table with 2 columns: Parameter (pH, Temperature, Background light, HOLD, Auto. switch off, Selection unit, Housing protection type, Batteries) and Specification.

Table with 2 columns: Action (Replacing the electrode, Storage of the electrode) and Description (rinsing, 3M KCl solution).

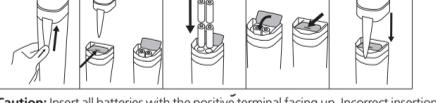Exchange batteries

Caution: Insert all batteries with the positive terminal facing up. Incorrect insertion of the batteries could result in damage to the meter.

Note: Make sure that the battery cover is completely closed to ensure water tightness.

Error messages and procedure

Table with 3 columns: Reason (Err 1, Err 2, Err 3) and Procedure (calibration sequence, old solution, key pressed too early).

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Zubehör

Table with 2 columns: Art.-Nr. (part number) and Beschreibung (description of accessories like buffer solution, tablets, beaker).

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Table with 3 columns: Calibration mode, Error code (Err 1, Err 2), Reason, and Procedure.

Note: In case of heavy contamination, carefully remove the sensor protection and rinse the electrode.

Accessories

Table with 2 columns: Art.-No. (part number) and Description (description of accessories).

Warranty

The warranty period for the SD10 pH is 2 years from the date of purchase and 6 months for the electrode.

Complete instructions with accessories list and further error messages can be found at www.lovbond.com (Download Centre).

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